

## JET BIOFIL® Super Clear Microcentrifuge Tubes



*Clearer Cheaper*

Biofil® Super clear tubes are made from an ultra-clear polypropylene that provides a clear view of the pellet. Polypropylene tubes feature a large cap to allow for easy opening and labeling. Flat caps and frosted side writing surface for easy labeling. Tight-fitting cap minimizes risk or popping OPEN due to heat.

**Avoid any possible improper manipulation, please read the instruction & note below carefully:**

- The maximum RCF is 16,000 g.
- Tubes are made by ultra super clear material. After heating samples, please wait the temperature down to RT then centrifuge.
- If centrifuge time is more than 15 minutes, please keep the centrifuging temperature in 4 °C.
- Don't over load samples more than 1.5 ml.
- Don't expose tubes in hood UV light too long.

*Please see more technical information  
on our website: [www. biotang.com](http://www.biotang.com)*

